

Installation Instructions Stands

KERN SXC, IXC

Type TSXC-A01-A
TIXC-A01-A
TIXC-A02-A

Version 1.0
2023-06
GB



TSXC-A01-A_TIXC-A01-A_TIXC-A02-A-IA-e-2310



Stands
Version 1.0 2023-06
Installation Instructions

Contents


1 Technical data..... 3
1.1 Dimensions 3
2 Scope of delivery 6
3 Installation..... 7

1 Technical data




KERN	TSXC-A01-A	TIXC-A01-A	TIXC-A02-A
Stand height (mm)	330	330	600

1.1 Dimensions




- Dimensions in mm
- Drawings on the following page

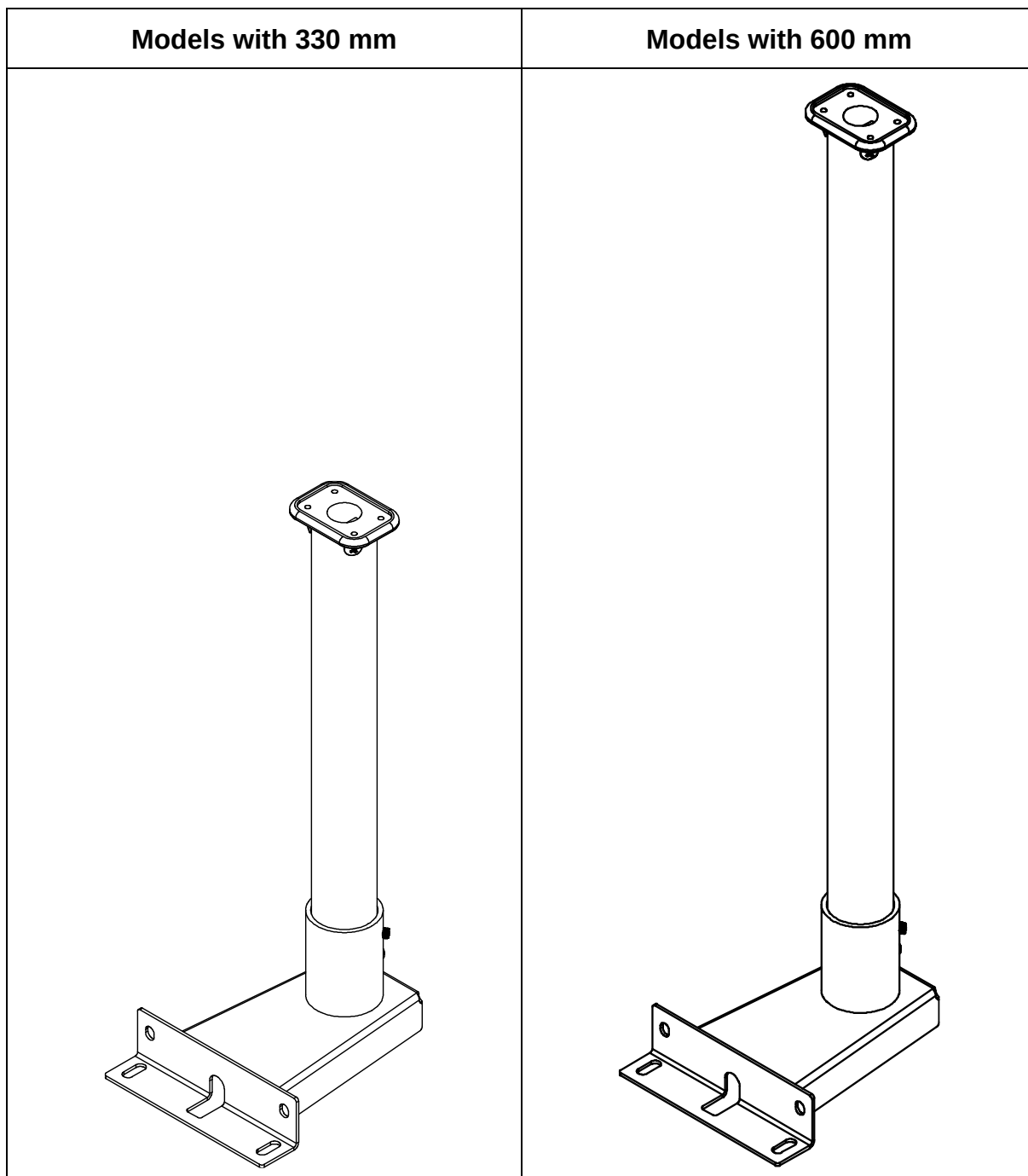







Models with 330 mm:

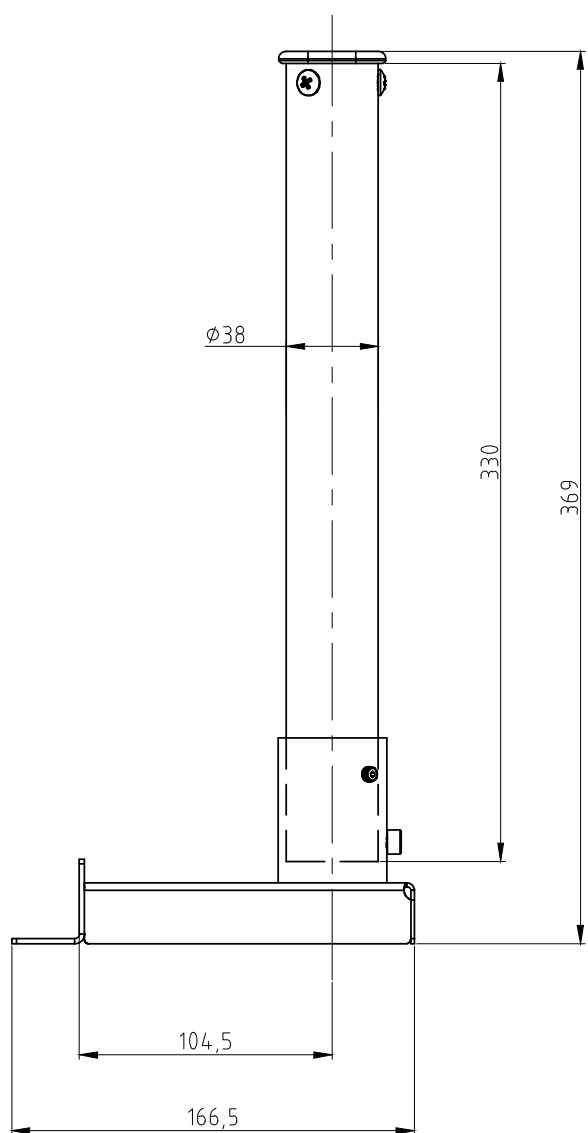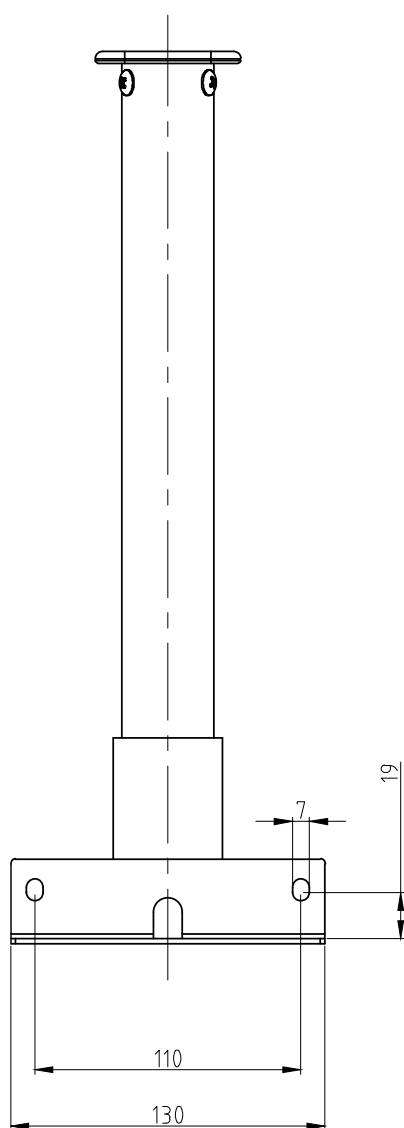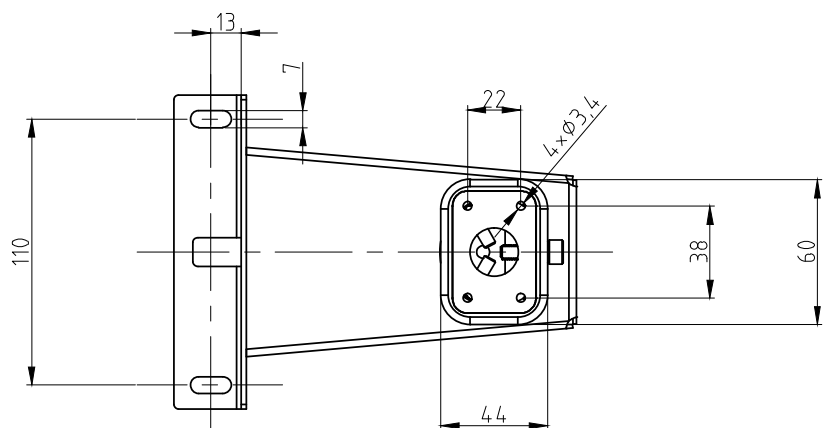

Models with 600 mm:

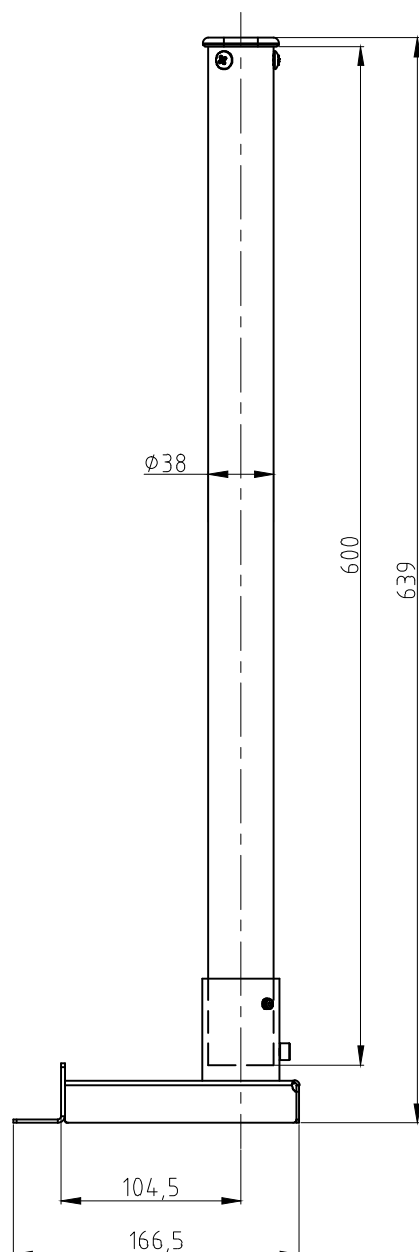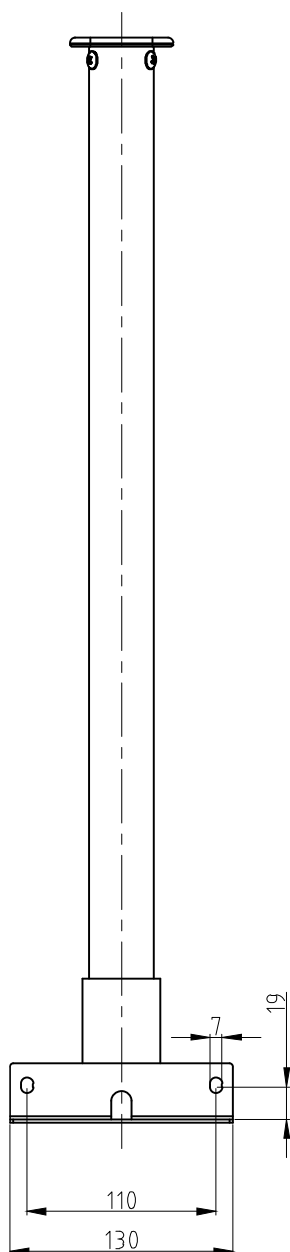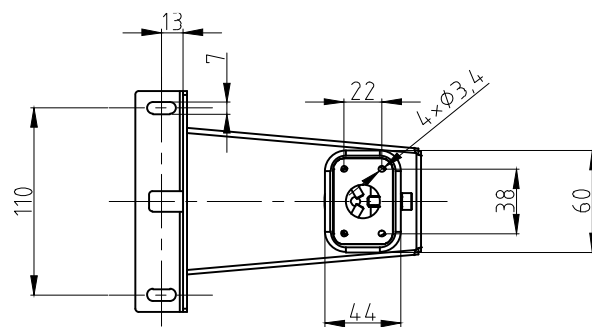

2 Scope of delivery

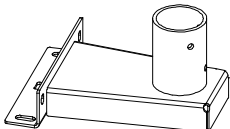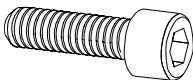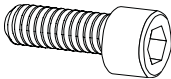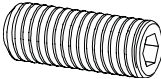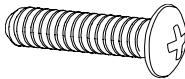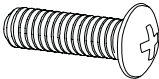
 A Stand foot 1 x	 B Stand tube 1 x	 C Support for display unit 1 x
 S1 M6 x 16 4 x	 S2 Spring washer 4 x	 S3 Washer 4 x
 S4 M6 x 12 1 x	 S5 M6 x 10 2 x	 S6 M5 x 16 3 x
 S7 M4 x 10 4 x		

3 Installation

Assembling the stand foot to the platform:


1. Place the stand foot (A) centrally over the perforations on the platform side.
2. First plug the spring washer (S2) and then the washer (S3) on the screw (S1) (4 x).
3. Screw down the stand foot.

Make sure that no cables are jammed or damaged.

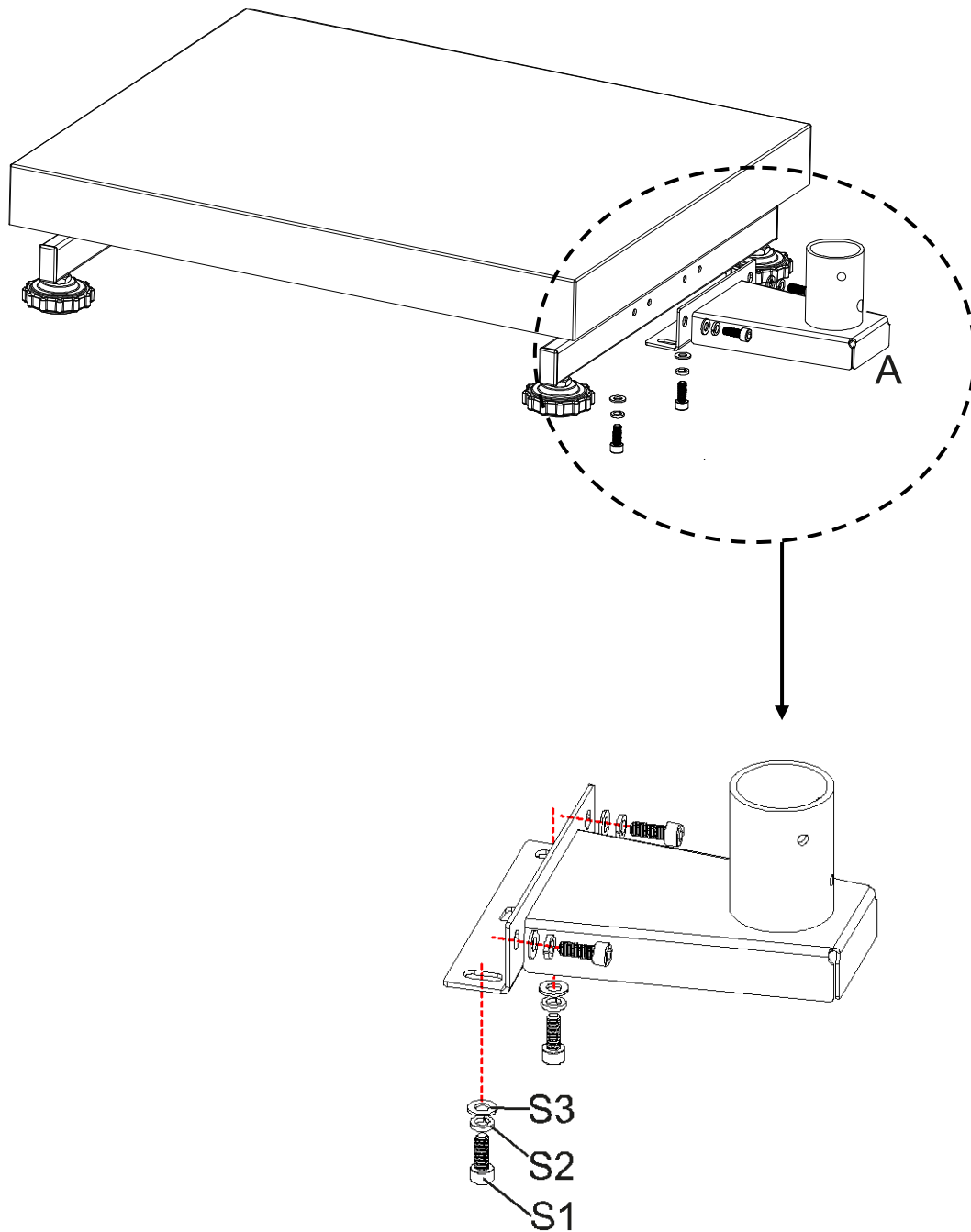

Assembling the stand tube to the stand foot:

4. Unscrew the screws **S4** and **S5** from the stand tube.
5. Push the stand tube **(B)** into the stand foot. Ensure that the perforation of the stand tube matches with the perforation **S4**.
6. Screw down again the screws **S5** and fasten the stand tube using screw **S4**.

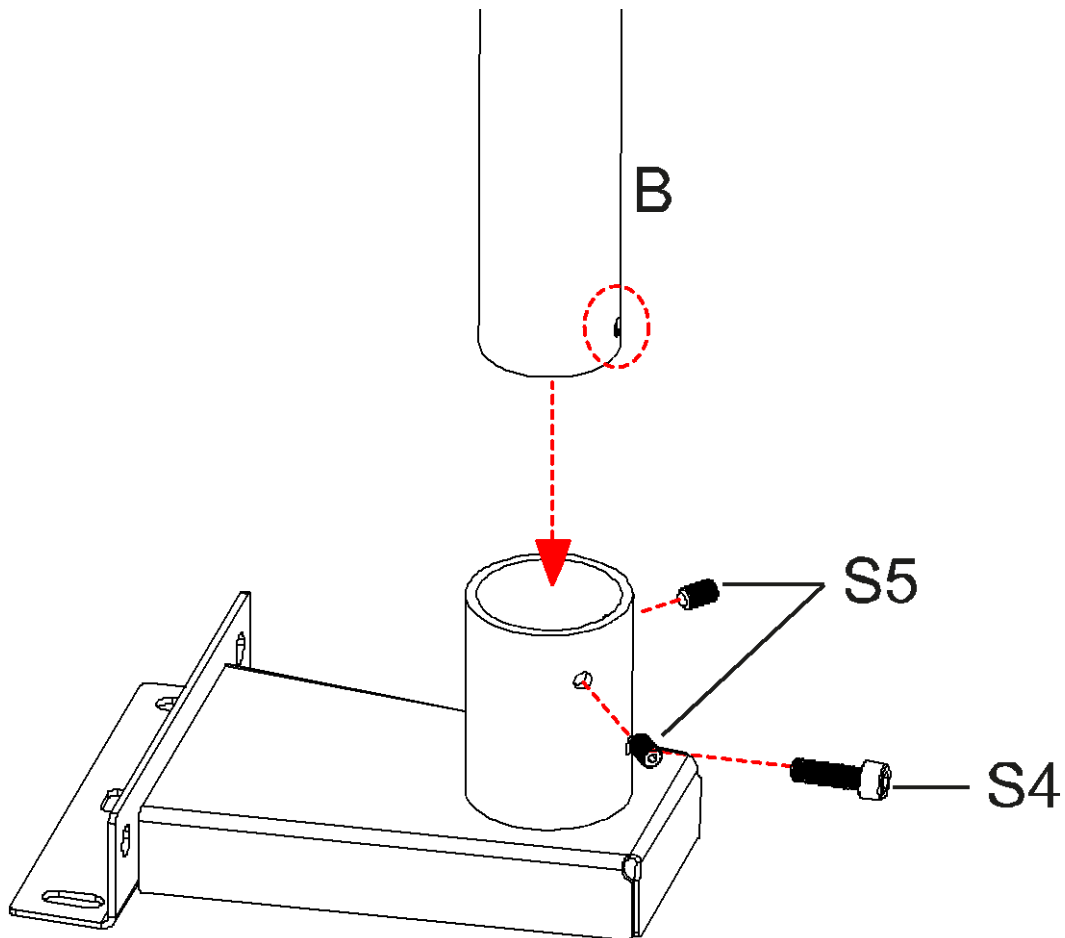

Assembling the display unit to the stand tube:

7. Push the support of the display unit (**C**) into the stand tube and fasten it using screws **S6**.
8. Fasten the display unit to the display unit's support using screws **S7**.

