

User instructions

Microscope camera

KERN ODC-85

ODC 851, ODC 852

Version 1.3
09/2023





KERN ODC-85

Version 1.3 09/2023

User instructions

Microscope camera

Before use

You should ensure that the device is not exposed to direct sunlight, temperatures which are too high or too low, vibrations, dust or a high level of humidity.

The ideal temperature range is between 0 and 40°C and a relative humidity of 85% should not be exceeded.

Always make sure that you use an approved power cable. Thus, possible damages by reason of the development of overheating (fire hazard) or an electric shock can be prevented.

Do not open the housing and touch the internal component. There is the risk of damaging them and affecting the functionality of the camera.

In order to carry out cleanings always disconnect the power cable from the camera.

Always keep the sensor clear from dust and do not touch it. Otherwise, there is the risk of affecting the microscopic image. In case of non-use always attach the protective covers.

Technical data

Model KERN	Resolution	Interface	Sensor	Frame rate	Colour / Monochrome	Supported operating systems
ODC 851	2 MP	HDMI, USB 2.0, SD	1/2,8" CMOS	30 - 60 fps	Colour	Win, XP, Vista, 7, 8, 10
ODC 852	5 MP	HDMI, USB 2.0, SD, WiFi	1/1,8" CMOS	25 - 60 fps	Colour	Win, XP, Vista, 7, 8, 10

Scope of delivery

- Microscope camera
- HDMI cable
- USB cable (ODC 851)
- SD card
- WiFi adapter (ODC 852)
- Object micrometre for calibration
- Software CD

Free download:

www.kern-sohn.com > DOWNLOADS > SOFTWARE > Microscope VIS Pro

- Eyepiece adapter (\varnothing 23,2 mm)
- Adjustment rings (\varnothing 30,0 mm + \varnothing 30,5 mm) for eyepiece adapter -
USB mouse
- Power supply

Mounting

1. Remove the black cover at the bottom of the camera.
2. The thread, where the cover was attached, is a standardised C-mount thread. Thus, there are special C mount adapters needed for the connection to a microscope.
3. For the mounting to the microscope the C mount adapter is attached to the connection point of the microscope. After that the camera must be screwed onto the C mount adapter.

Important:

The choosing of the right C mount adapter depends on the used microscope model. It has to be an adapter, which is adjusted to the construction of the microscope and recommended by the manufacturer as appropriate for the relevant microscope.

4. If necessary, adjust the microscope according to the trinocular usage (with help of the trino toggle rod / trino toggle wheel).

The KERN ODC-85 series ensures to carry out digital microscopy either with help of an HDMI connection directly on the screen or with help of an USB 2.0/WiFi connection (via software).

Screen connection (HDMI)

1. Establish HDMI connection via HDMI cable and switch on the camera by the power button.
2. Stick SD card into SD port on the camera.
3. Connect USB mouse to USB port of the camera.
4. As soon as the image transmission has been launched, the cursor appears on the screen. When moved to the edge of the screen, it folds out some editing menus and further control elements (e. g. for data storage).
5. Short function explanations are integrated to each of the selectable control elements (in English).

PC connection ODC 851 (USB 2.0)

1. Establish USB connection via USB cable and switch on the camera by the power button.
2. Installing the software with help of the CD / download.
3. The software-internal "User Guide" includes all information and instructions about the operation of the software or of digital microscopy.

PC connection ODC 851 (WiFi)

1. Plug in the WiFi adapter to the USB port of the camera and switch on the camera by the power button.
2. With activated WiFi antenna of the PC, the hotspot of the camera is shown in the network settings:

„XFCAM1080PHB_#“ Key: 12345678

3. Installing the software with help of the CD / download.
4. The software-internal "User Guide" includes all information and instructions about the operation of the software or of digital microscopy.