

Installation Instructions Stands

KERN SXC, IXC

Type TSXC-A01-A
TIXC-A01-A
TIXC-A02-A

Version 1.0
2023-06
GB



TSXC-A01-A_TIXC-A01-A_TIXC-A02-A-IA-e-2310



Stands

Version 1.0 2023-06

Installation Instructions

Contents


1	Technical data	3
1.1	Dimensions	3
2	Scope of delivery	6
3	Installation	7

1 Technical data




KERN	TSXC-A01-A	TIXC-A01-A	TIXC-A02-A
Stand height (mm)	330	330	600

1.1 Dimensions




- Dimensions in mm
- Drawings on the following page

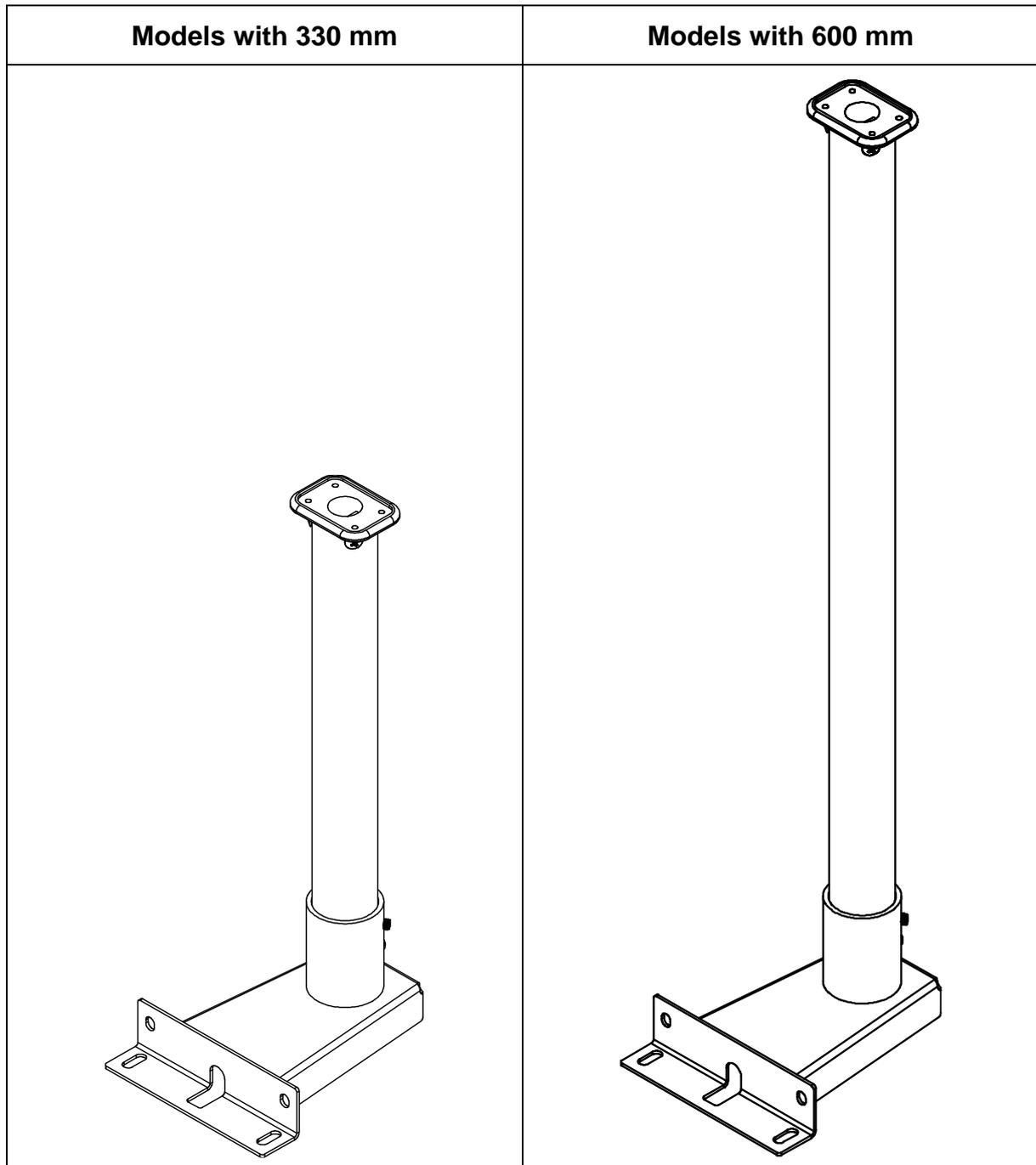







Models with 330 mm:

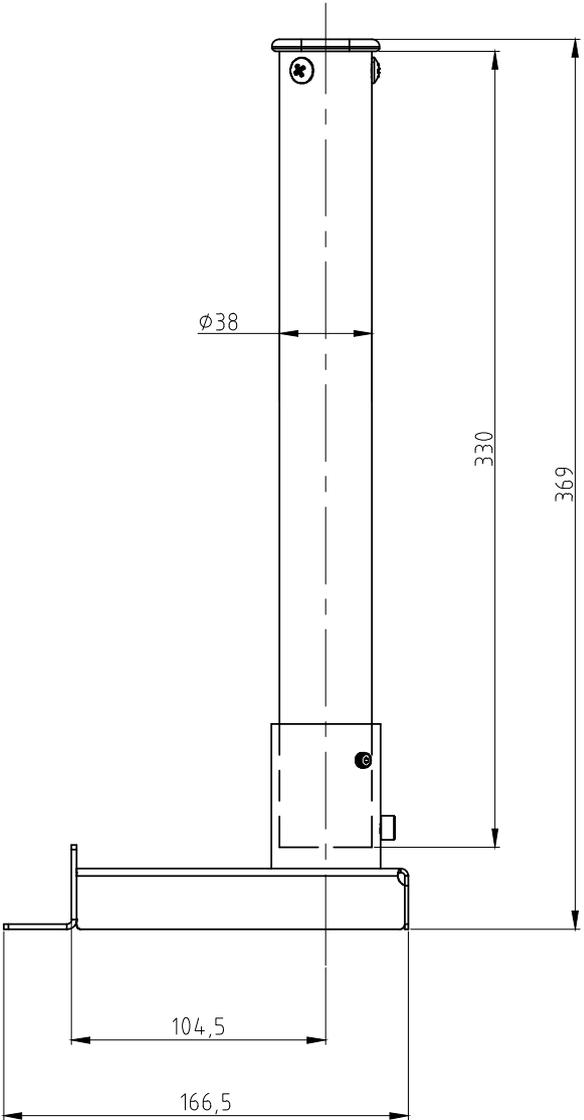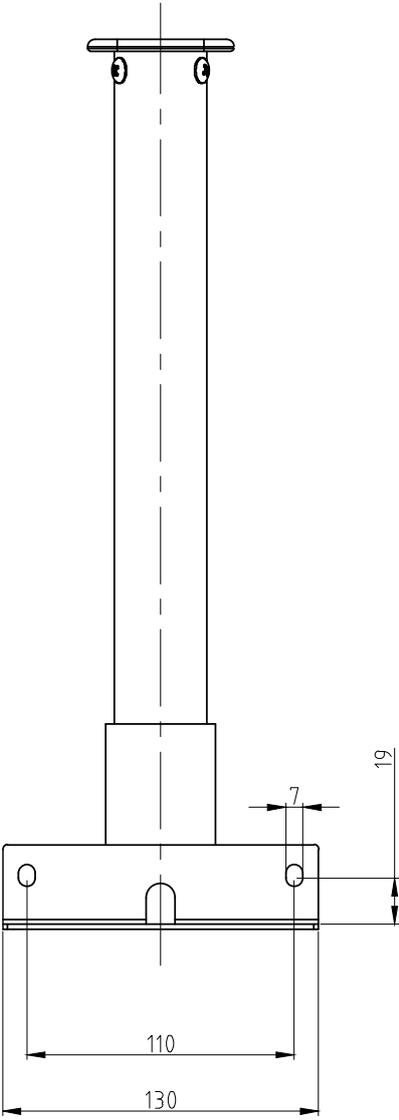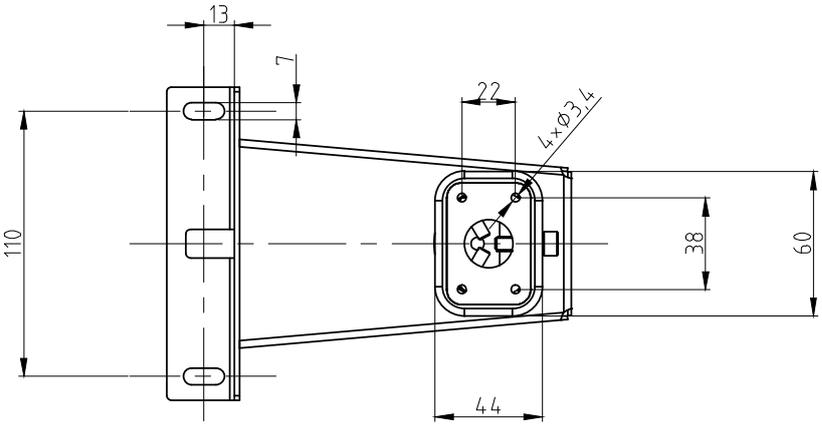

Models with 600 mm:

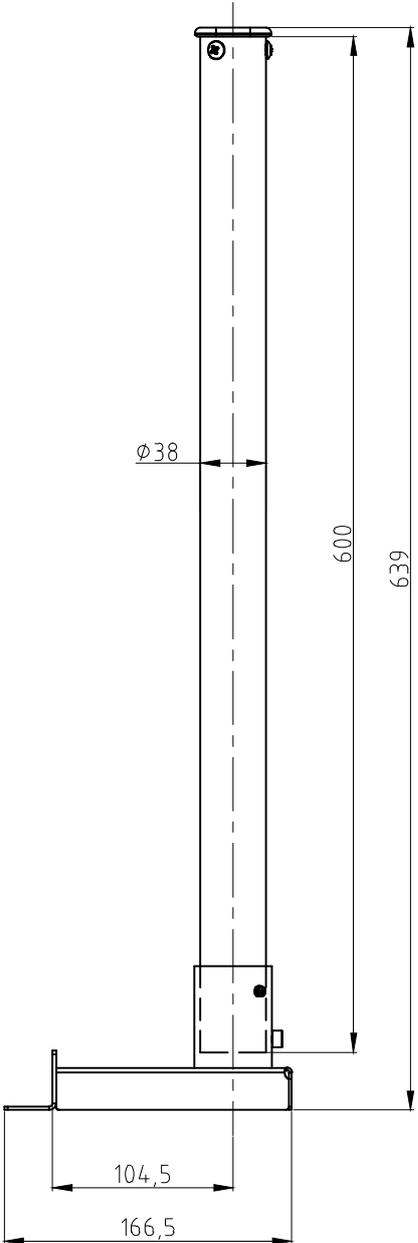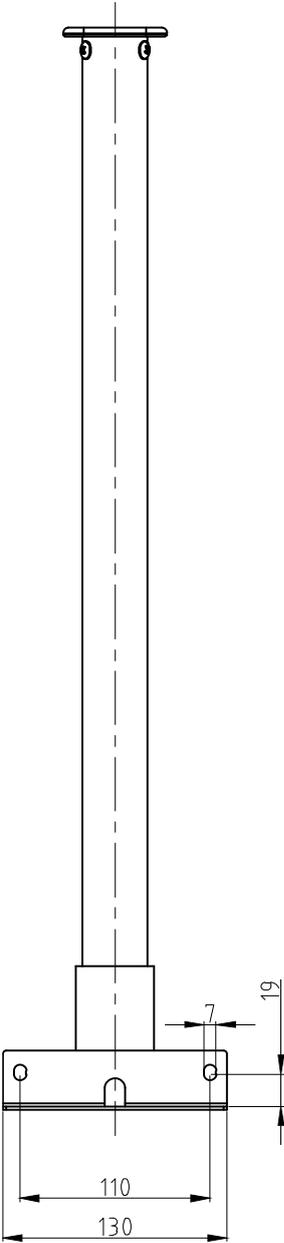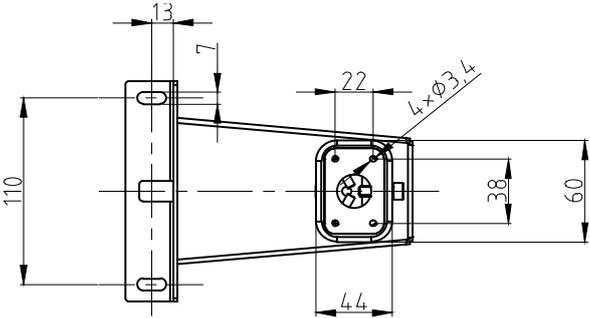

2 Scope of delivery

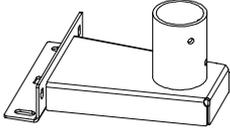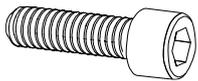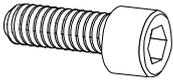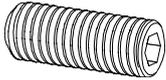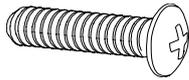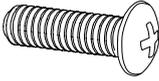
 <p>A</p> <p>Stand foot</p> <p>1 x</p>	 <p>B</p> <p>Stand tube</p> <p>1 x</p>	 <p>C</p> <p>Support for display unit</p> <p>1 x</p>
 <p>S1</p> <p>M6 x 16</p> <p>4 x</p>	 <p>S2</p> <p>Spring washer</p> <p>4 x</p>	 <p>S3</p> <p>Washer</p> <p>4 x</p>
 <p>S4</p> <p>M6 x 12</p> <p>1 x</p>	 <p>S5</p> <p>M6 x 10</p> <p>2 x</p>	 <p>S6</p> <p>M5 x 16</p> <p>3 x</p>
 <p>S7</p> <p>M4 x 10</p> <p>4 x</p>		

3 Installation

Assembling the stand foot to the platform:


1. Place the stand foot (**A**) centrally over the perforations on the platform side.
2. First plug the spring washer (**S2**) and then the washer (**S3**) on the screw (**S1**) (4 x).
3. Screw down the stand foot.

Make sure that no cables are jammed or damaged.

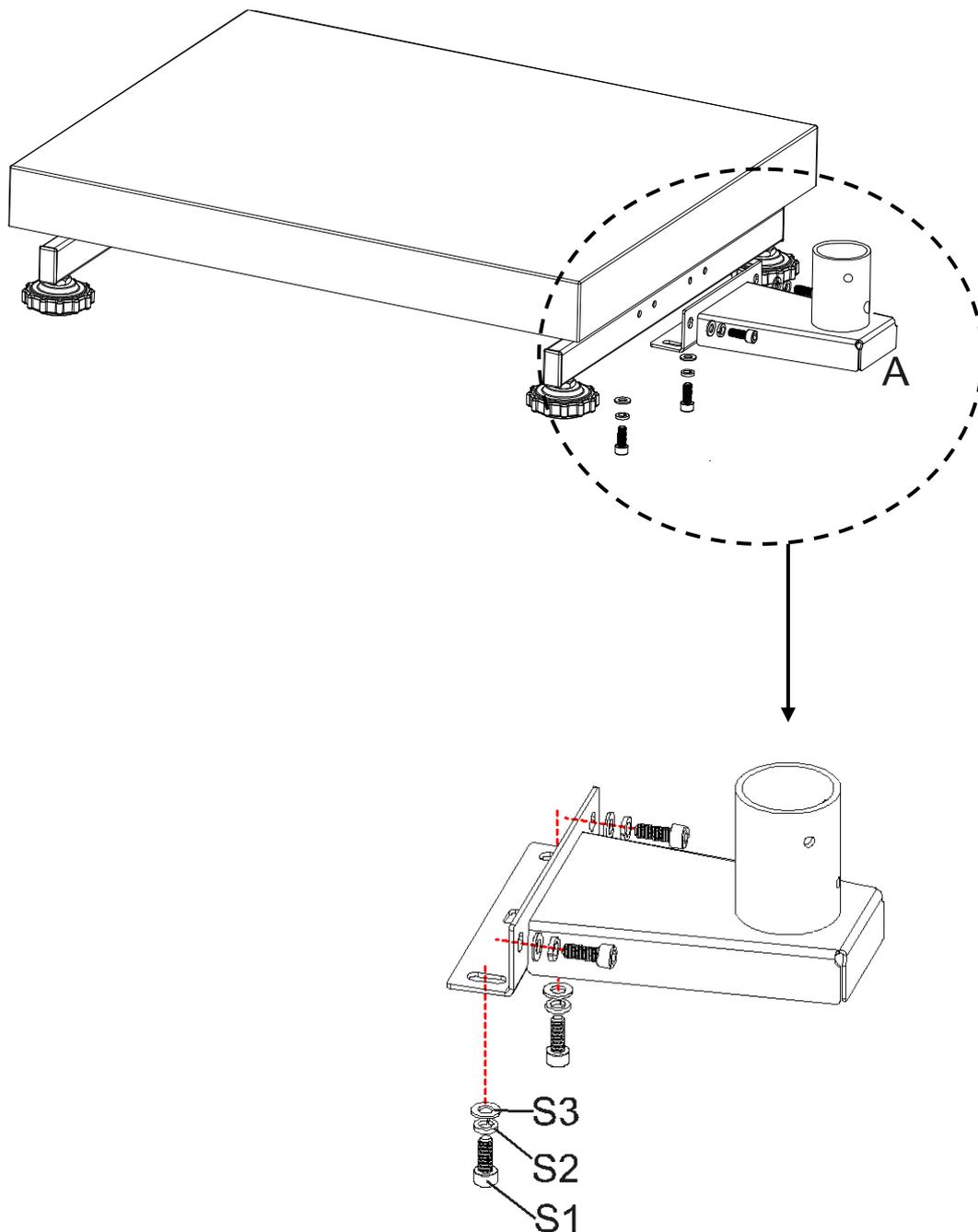

Assembling the stand tube to the stand foot:

4. Unscrew the screws **S4** and **S5** from the stand tube.
5. Push the stand tube (**B**) into the stand foot. Ensure that the perforation of the stand tube matches with the perforation **S4**.
6. Screw down again the screws **S5** and fasten the stand tube using screw **S4**.

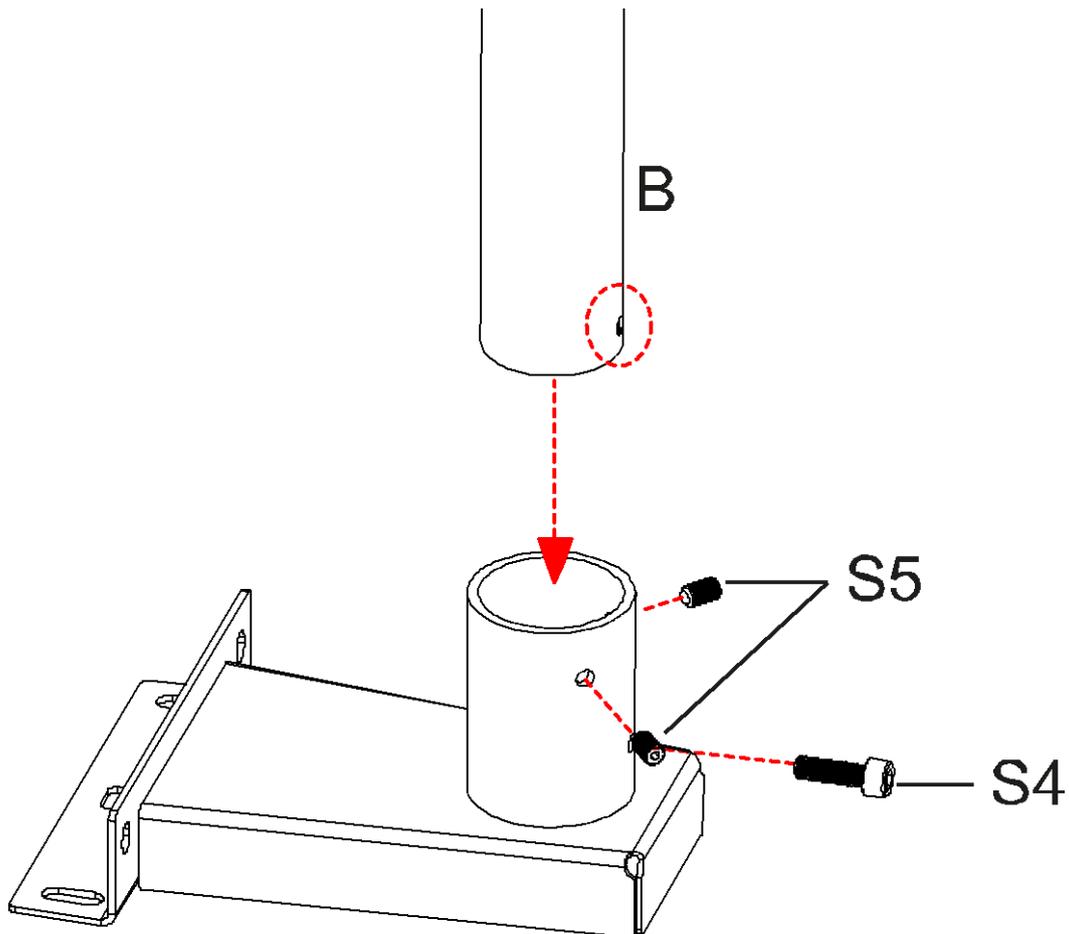

Assembling the display unit to the stand tube:

7. Push the support of the display unit (**C**) into the stand tube and fasten it using screws **S6**.
8. Fasten the display unit to the display unit's support using screws **S7**.

