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Operating Instructions

1 Insert handle – except when the scale is hung up for weighing.

2 Adjust zero position – with the scale in measuring position.

3 Tare: Attach empty container or other tare if required and adjust zero position.

4 Position drag pointer to zero – if not in use simply remove. The application of the drag pointer might very slightly influence measurement precision.

5 Neither load scale abruptly nor in excess of 130 % of capacity. Burst tests limited to 30 kg/300 N (possible damage by setback). This product is not to be used by children.

Options

6 Exchange of lower suspension (accessories): see special instructions overleaf.

7 Transformation to pressure dynamometer with accessory 8.004.

Maintenance

No particles should get into the scale. Wash away salt water and dirt. Otherwise the scale needs no maintenance, is non corrosive and has an unlimited life with constant precision.

Precision

Maximum tolerance ± 0.3 % of load (zero position exactly adjusted, temperature 10 – 30 °C/ 50 – 86 °F and using neither drag pointer nor pressure-set).

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English

Exchange of lower suspension of Macro Scales with optional accessories

ATTENTION: Should this instruction not be observed, damages may occur which will not be covered by warranty.

1 Keep on turning the adjustment screw to the left (A) until the inner plastic tube falls down to its lower stop.

2 With the scale upside-down, hold the inner plastic tube sticking out of the outer aluminum tube with your hand (A). Unscrew the hexagonal nut of the hook with a spanner wrench (12 mm) by one turn (B). (ATTENTION: Do not counter-hold the outer aluminum tube when unscrewing the hexagonal nut with the wrench. The torsion force might damage the small marker in the slit of the scale.) Unscrew the hexagonal nut of the hook completely by hand. (ATTENTION: Should the hexagonal nut idle and not allow unscrewing or tightening, an inner cam in the plastic tube might be cut off and needs service by the PESOLA factory.)

3 Slightly press the inner plastic tube into the outer aluminum tube against the force of the inner spring (A) and rotate the threaded bolt M8 until it clicks into its place (B), showing its maximum length. Existing discs and/or washers shall be used again.

4 Maintain a minimum pressure on the inner plastic tube (A) so as to keep the threaded bolt M8 in its firm position. (Optional: Possibly apply some weak glue on the threaded bolt M8.) Fully screw the hexagonal nut M8 of the optional lower suspension (adaptation set 8.018, heavy duty clamp 8.010, replacement hook 8.128, cap nut M8 of pressure set 8.004 or other) on the threaded bolt by hand only (B).

5 Fully extend the inner plastic tube and counter-hold it with your hand (A). (ATTENTION: Do not counter-hold the outer aluminum tube when tightening the hexagonal nut M8. Torsion forces may damage the small marker in the slit.) Smoothly tighten the hexagonal nut M8 with a 12 mm spanner wrench without using excessive force (B).

6 Still holding the scale upside down, twist the adjustment screw to the right (A). The inner tube will slowly move back into the outer aluminum tube. For your next weighing job precisely adjust the zero point with the scale in measuring position.

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