PX Series Balances User Guide

TABLE OF CONTENT

EN	
1. INTRODUCTION	3
2. SAFETY INFORMATION	3
3. INSTALLATION	4
4. OPERATION	5
5. MAINTENANCE	6
6. TECHNICAL DATA	6
ES	
1. INTRODUCCIÓN	8
2. INFORMACIÓN SOBRE SEGURIDAD.....	8
3. INSTALLATION	9
4. FUNCIONAMIENTO	10
5. MANTENIMIENTO	11
6. DATOS TÉCNICOS.....	11
FR	
1. INTRODUCTION	13
2. INFORMATIONS DE SECURITE.....	13
3. INSTALLATION	14
4. OPERATION	15
5. ENTRETIEN	16
6. DONNÉES TECHNIQUES	16
DE	
1. EINFÜHRUNG	18
2. SICHERHEITSINFORMATIONEN	18
3. INSTALLATION	19
4. BETRIEB	20
5. WARTUNG	21
6. TECHNISCHE DATEN	22
IT	
1. INTRODUZIONE	23
2. INFORMAZIONI SULLA SICUREZZA	23
3. INSTALLAZIONZ	24
4. FUNZIONAMENTO	25
5. MANUTENZIONE	26
6. DATI TECNICI	27
PL	
1. WPROWADZENIE	28
2. INFORMACJE DOTYCZĄCE BEZPIECZEŃSTWA	28
3. INSTALACJA	29
4. DZIAŁANIE	30
5. KONSERWACJA	31
6. DANE TECHNICZNE	32
SV	
1. INLEDNING	33
2. SÄKERHETSFÖRESKRIFTER.....	33
3. INSTALLATION	34
4. DRIFT	35
5. UNDERHÅLL	36
6. TEKNISKA DATA	36

NL	
1. INLEIDING.....	38
2. VEILIGHEIDSGEGEVENS.....	38
3. INSTALLATIE	39
4. GEBRUIK	40
5. ONDERHOUD	41
6. TECHNISCHE GEGEVENS	42
DA	
1. INDLEDNING	43
2. OPLYSNINGER VEDRØRENDE SIKKERHED.....	43
3. INSTALLATION	44
4. BETJENING	45
5. VEDLIGEHOLDELSE	46
6. TEKNISKE DATA.....	46
PT	
1. INTRODUÇÃO	48
2. INFORMAÇÕES DE SEGURANÇA.....	48
3. INSTALAÇÃO	49
4. FUNCIONAMENTO.....	50
5. MANUTENÇÃO.....	51
6. DADOS TÉCNICOS	52
CZ	
1. ÚVOD	53
2. BEZPEČNOSTNÍ INFORMACE.....	53
3. INSTALACE	54
4. PROVOZ.....	55
5. ÚDRŽBA.....	56
6. TECHNICKÉ ÚDAJE.....	56
HU	
1. BEVEZETÉS	58
2. BIZTONSÁGI TUDNIVALÓK.....	58
3. ÜZEMBE HELYEZÉS.....	59
4. HASZNÁLAT.....	60
5. KARBANTARTÁS.....	61
6. TECHNICAL DATA.....	61

1. INTRODUCTION

Thank you for purchasing our product. This user guide will help you install your balance and use it safely. Please read it carefully before installation and operation. For detailed instructions, please visit www.ohaus.com/products.

2. SAFETY INFORMATION

This manual contains installation, operation and maintenance instructions for PX Series balance. Please read this manual completely before installation and operation.

Definition of Signal Warnings and Symbols

Safety notes are marked with signal words and warning symbols. These show safety issues and warnings. Ignoring the safety notes may lead to personal injury, damage to the instrument, malfunctions and false results.

WARNING For a hazardous situation with medium risk, possibly resulting in injuries or death if not avoided.

CAUTION For a hazardous situation with low risk, resulting in damage to the device or the property or in loss of data, or injuries if not avoided.

Attention Note For important information about the product

Note For useful information about the product

Warning Symbols

General Hazard

Electrical Shock Hazard

Alternating Current

Direct Current

Safety Precautions

CAUTION: Read all safety warnings before installing, making connections, or servicing this equipment. Failure to comply with these warnings could result in personal injury and/or property damage. Retain all instructions for future reference.

- Verify that the AC adapter's input voltage range and plug type are compatible with the local AC main power supply.
- Make sure that the power cord does not pose a potential obstacle or tripping hazard.
- Do not position the balance such that it is difficult to reach the power connection.
- The balance is for indoor use only. Do not operate the equipment in hazardous or unstable environments.
- Operate the equipment only under ambient conditions specified in these instructions.
- Do not drop loads on the pan.
- Use the balance only in dry locations.
- Disconnect the equipment from the power supply when cleaning.
- Use only approved accessories and peripherals.
- Service should only be performed by authorized personnel.

Intended Use

Use the instrument exclusively for weighing as described in the operating instructions.

Any other type of use and operation beyond the limits of technical specifications without written consent from OHAUS, is considered as not intended.

This instrument complies with current industry standards and the recognized safety regulations; however, it can constitute a hazard in use.

If the instrument is not used according to these operating instructions, the intended protection of the instrument may be compromised and OHAUS assumes no liability.

3. INSTALLATION

3.1 Select the Location

Avoid heat sources

Rapid temperature changes

Air current or excessive vibrations

Allow sufficient space

3.2 Leveling

Be sure the balance is level before it is used or after its location is changed.

The PX balance has a level bubble in a small round window beside the display.

To level the balance, adjust the 4 leveling feet until the bubble is centered in the circle.

Please refer to the right figure for leveling.

3.3 Connecting Power and Acclimatising the Balance

Connect the DC output connector to the power receptacle on the rear of the balance. Then connect the AC adapter plug to a suitable electrical outlet.

Acclimatising

It is suggested that the balance should not be used until it has been connected to power and acclimatised to the environment for a certain period of time. In the case of a balance with the precision above 0.1 mg, the acclimatisation time should be 1.5 hours; in the case of balance with the precision of 0.01 mg, the acclimatisation time should be more than 4 hours.

3.4 Connecting the Interface

The PX balance has two data interfaces, RS232 and USB.

Use the RS-232 port to connect either to a computer or a printer with a standard (straight-through) serial cable. Use the USB port to connect to a computer with a USB 2.0 Type A to Type B cable.

Interface connections on the rear of the balance

USB

RS232

USB: Used to connect to PC only

RS232: Used to connect to PC or Printer

Note: For Connecting, Configuring and Testing the Printer / Computer Interface, see the Printing section.

3.5 Initial Calibration

For best results, the instrument's calibration should be checked prior to first use. If adjustment is needed, refer to the Calibration section of the instruction manual.

4. OPERATION





4.1 Overview of Controls and Display

The PX balance has a 2-line backlit display.

CONTROLS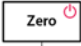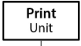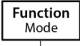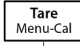

CONTROL FUNCTIONS

Button	Zero 	Print Unit 	Function Mode 	Tare Menu-Cal 
Primary Function (Short Press) 	On / Zero <ul style="list-style-type: none"> If the balance is Off, turns on the balance. If the balance is On, sets zero. 	Print <ul style="list-style-type: none"> Sends the current displayed value to the serial interface. 	Function <ul style="list-style-type: none"> Operation is dependent on the application mode. 	Tare <ul style="list-style-type: none"> Performs tare operation.
Secondary Function (Press and Hold)  (Extended Press)	Off <ul style="list-style-type: none"> Zeros the current value. None 	Unit <ul style="list-style-type: none"> Changes weighing units. None 	Mode <ul style="list-style-type: none"> Changes application mode. None 	Menu-Cal <ul style="list-style-type: none"> Enters the main menu. Calibration is the first sub-menu. Views the preset Tare value.
Menu Function (Short Press) 	Yes <ul style="list-style-type: none"> Accepts the current (blinking) setting on the display. 	No <ul style="list-style-type: none"> Rejects the current (blinking) setting on the display. Increases a value being entered. 	Back <ul style="list-style-type: none"> Reverts back to previous menu item. Decrements a value being entered. 	Exit <ul style="list-style-type: none"> Immediately exits the sub-menu. Aborts a calibration in progress.

DISPLAY

Net (NET)
 Stability (*)
 Negative (-)
 Centre of Zero (>0<)
 Preset tare (PT)
 Instructional Messages

Result Field:
 Information varies by application
 Unit

4.2 Principal Functions

Zeroing:	Press Zero to zero the balance.
Taring:	With no load on the pan, press Zero to set the display to zero. Place an empty container on the pan and press Tare . Add material to the container and its net weight is displayed. After the container and the objects are removed, the tare value will be displayed as a negative number. Press Tare to clear.
Weighing:	First press Zero to set the display to zero. Place an object on the pan. The display indicates the gross weight.
Dot-matrix Display:	The relevant data in the specific application mode are shown in the dot-matrix display area.

5. MAINTENANCE

5.1 Cleaning

WARNING: Disconnect the balance from the power supply before cleaning. Make sure that no liquid enters the interior of the balance.

Clean the balance at regular intervals.

Housing surfaces may be cleaned with a lint-free cloth slightly dampened with water or a mild cleaning agent.

Glass surfaces may be cleaned with a commercial glass cleaner.

Attention: Do not use solvents, harsh chemicals, ammonia or abrasive cleaning agents.

5.2 Service Information

For troubleshooting common problems, download the user manual from www.Ohaus.com. If the issue persists, please visit our website for available support and service centers at <https://www.ohaus.com/en-us/support/service-centers>.

6. TECHNICAL DATA

The technical data is valid under the following conditions:

- Indoor use only
- Altitude: Up to 2000 m
- Operating temperature: 5°C to 40°C
- Operating temperature for optimum weighing performance: 10°C to 30°C
- Humidity: maximum relative humidity 80% for temperatures up to 30°C, decreasing linearly to 50% relative humidity at 40°C
- Electrical supply: 12VDC, 0.5A. (For use with certified or approved power supply, which must have a SELV and limited energy output.)
- Mains supply voltage fluctuations: up to $\pm 10\%$ of the nominal voltage
- Installation category II
- Pollution degree: 2

Detailed product specifications can be found in the user manual and can be downloaded from www.Ohaus.com/products.

LIMITED WARRANTY

Ohaus products are warranted against defects in materials and workmanship from the date of delivery through the duration of the warranty period. During the warranty period Ohaus will repair, or replace any component(s) that proves to be defective at no charge, provided that the product is returned, freight prepaid, to Ohaus.

This warranty does not apply if the product has been damaged by accident or misuse, exposed to radioactive or corrosive materials, has foreign material penetrating to the inside of the product, or as a result of service or modification by other than Ohaus. In lieu of a properly returned warranty registration card, the warranty period shall begin on the date of shipment to the authorized dealer. No other express or implied warranty is given by Ohaus Corporation. Ohaus Corporation shall not be liable for any consequential damages.

As warranty legislation differs from state to state and country to country, please contact Ohaus or your local Ohaus dealer for further details.

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