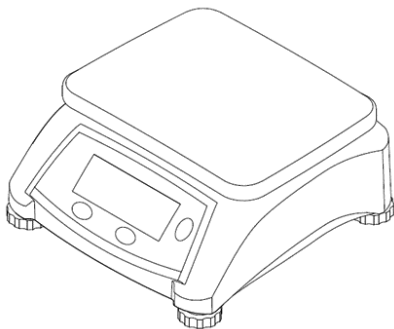
Valor™ 1000 Series User Guide

REVISION HISTORY

Date	Version	Description
February 14 2023	D	We incorporate the product specification table into the instruction manual in an effort to reduce printed materials. Please download the user instruction manual at ohaus.com/products or use your smartphone to scan the QR code on the front page of this manual.

TABLE OF CONTENTS

EN	
1.	INTRODUCTION 3
2.	SAFETY INFORMATION 3
3.	INSTALLATION 4
4.	OPERATION 5
5.	MAINTENANCE 6
6.	TECHNICAL DATA 7
ES	
1.	INTRODUCCIÓN 8
2.	INFORMACIÓN DE SEGURIDAD 8
3.	INSTALACIÓN 9
4.	OPERACIÓN 10
5.	MANTENIMIENTO 11
6.	DATOS TÉCNICOS 12
FR	
1.	INTRODUCTION 13
2.	INFORMATIONS DE SECURITE 13
3.	INSTALLATION 14
4.	FONCTIONNEMENT 15
5.	ENTRETIEN 16
6.	DONNÉES TECHNIQUES 17
DE	
1.	EINLEITUNG 18
2.	SICHERHEITSINFORMATIONEN 18
3.	EINRICHTUNG 19
4.	BEDIENUNG 20
5.	WARTUNG 21
6.	TECHNISCHE DATEN 22
IT	
1.	INTRODUZIONE 23
2.	INFORMAZIONI DI SICUREZZA 23
3.	INSTALLAZIONE 24
4.	FUNZIONAMENTO 25
5.	MANUTENZIONE 27
6.	DATI TECNICI 27
PT	
1.	INTRODUÇÃO 28
2.	INFORMAÇÕES DE SEGURANÇA 28
3.	INSTALAÇÃO 29
4.	FUNIONAMENTO 30
5.	MANUTENÇÃO 31
6.	ESPECIFICAÇÕES TÉCNICAS 32

SV	
1.	INLEDNING..... 33
2.	SÄKERHETSINFORMATION..... 33
3.	INSTALLATION..... 34
4.	DRIFT..... 35
5.	UNDERHÅLL..... 36
6.	TEKNISK DATA 37
NL	
1.	INLEIDING 38
2.	VEILIGHEIDSINFORMATIE 38
3.	INSTALLATIE 39
4.	WERKING 40
5.	ONDERHOUD 41
6.	TECHNISCHE GEGEVENS 42
DA	
1.	INDLEDNING 43
2.	SIKKERHEDSINFORMATION 43
3.	INSTALLATION..... 44
4.	BRUGSVEJLEDNING 45
5.	VEDLIGEHOOLD 46
6.	TEKNISK DATA 47
PL	
1.	WPROWADZENIE 48
2.	INFORMACJE DOTYCZĄCE BEZPIECZEŃSTWA 48
3.	INSTALACJA..... 49
4.	OBSŁUGA..... 50
5.	KONSERWACJA..... 51
6.	DANE TECHNICZNE 52
CS	
1.	ÚVOD 53
2.	BEZPEČNOSTNÍ INFORMACE 53
3.	INSTALACE 54
4.	OBSLUHA 55
5.	ÚDRŽBA..... 56
6.	TECHNICKÉ ÚDAJE 57
HU	
1.	BEVEZETÉS 58
2.	BIZTONSÁGI TUDNIVALÓK..... 58
3.	BESZERELÉS 59
4.	KEZELÉS 60
5.	KARBANTARTÁS..... 61
6.	MŰSZAKI ADATOK..... 62

1. INTRODUCTION

Thank you for purchasing our product. This user guide will help you install your scale and use it safely. Please read it carefully before installation and operation. For detailed instructions, please visit www.ohaus.com/products.

2. SAFTY INFORMATION

Definition of Signal Warnings and Symbols

Safety notes are marked with signal words and warning symbols. These show safety issues and warnings. Ignoring the safety notes may lead to personal injury, damage to the instrument, malfunctions and false results.

WARNING	For a hazardous situation with medium risk, possibly resulting in severe injuries or death if not avoided.
CAUTION	For a hazardous situation with low risk, resulting in damage to the device or the property or in loss of data, or minor or medium injuries if not avoided.
ATTENTION	For important information about the product. May lead to equipment damage if not avoided.
NOTE	For useful information about the product.

Warning Symbols

General hazard

Explosion hazard

Electrical shock hazard

Safety Precautions

CAUTION: Read all safety warnings before installing, making connections, or servicing this equipment. Failure to comply with these warnings could result in personal injury and/or property damage. Retain all instructions for future reference.

- Before connecting power, verify that the AC adapter's input voltage range and plug type are compatible with the local AC mains power supply.
- Do not position the equipment such that it is difficult to reach the power connection.
- Make sure that the power cord does not pose a potential obstacle or tripping hazard.
- Operate the equipment only under ambient conditions specified in these instructions.
- The equipment is for indoor use only.
- Do not operate the equipment in wet, hazardous or unstable environments.
- Do not allow liquids to enter the equipment.
- Do not load the equipment above its rated capacity.
- Do not drop loads on the platform.
- Do not place the equipment upside down on the platform.
- Use only approved accessories and peripherals.
- Disconnect the equipment from the power supply when cleaning.
- Service should only be performed by authorized personnel.

WARNING: Never work in an environment subject to explosion hazards! The housing of the instrument is not gas tight. (explosion hazard due to spark formation, corrosion caused by the ingress of gases).

CAUTION: Battery is to be replaced only by an authorized Ohaus service dealer. Risk of explosion can occur if the rechargeable battery is replaced with the wrong type or if it is not properly connected. Dispose of the rechargeable battery according to local laws and regulations.

Intended Use

This instrument is intended for use in businesses and light industry. It shall only be used for measuring the parameters described in these operating instructions. Any other type of use and operation beyond the limits of technical specifications, without written consent from OHAUS, is considered as not intended. This instrument complies with current industry standards and the recognized safety regulations; however, it can constitute a hazard in use. If the instrument is not used according to these operating instructions, the intended protection provided by the instrument may be impaired

3. INSTALLATION

3.1 Unpacking

Carefully remove your Valor 1000 scale and each of its components from the package. Save the packaging to ensure safe storage and transport. Please read the manual completely before installing and using the Valor 1000 scale to avoid incorrect operation.

Components included:

- Scale
- User Guide
- Stainless steel pan
- 4 dry batteries

3.2 Selecting the location

The location must be sturdy, flat and level. Avoid locations with excessive air current, vibration, heat sources or rapid temperature changes. Allow sufficient space around the instrument.

3.3 Power

Connect the AC Adaptor to the AC mains supply. Connect the plug to the DC jack on the bottom of the scale. The scale may be operated on the AC Adapter connected to mains supply or 4 alkaline batteries.

Note: The AC adapter maybe optional or included with the scale, depending on the model.

3.4 Rechargeable Battery Power

Allow the battery to charge for 12 hours before using the scale on battery power. The instrument can be operated during charging, and the battery is protected against overcharging. For maximum operating time, the battery should be charged at room temperature.

Note: Availability is depending on region as accessory.

WARNING: Only charge the battery when the ambient temperature is between 0 and 140 °F / 0 and 40 °C.

3.5 Levelling the Instrument

To level the instrument, adjust the feet/leveling wheel so the bubble is centered in the circle of the level indicator. Be sure the equipment is level each time its location is changed.

3.6 Initial Calibration

For best results, the instrument's calibration should be checked prior to first use. If adjustment is needed, refer to the Calibration section of the instruction manual.

4. OPERATION

Button	Functions
	On/Off – Press (when off): Turn the scale on – Long Press (when on): Turn the scale off Zero – Short Press (when on): Zero the scale Yes – Press (in Menu): Confirm (Enter)
	Tare – Short Press: Enter / clear a Tare value Menu – Long Press: Enter User Menu No – Press (in Menu): Move to next menu or selection

Item	Description
1	Battery charge symbol
2	Center of Zero symbol
3	Negative symbol
4	Stable weight symbol
5	NET symbol

Notes:

Short Press: Press less than 3 seconds.

Long Press: Press and hold for more than 3 seconds.

4.1 Using the Instrument

Weighing application:

1. If required, place an empty container on the pan and press Tare.
2. Add sample to the pan or container. The display shows the weight of the sample.

5. MAINTENANCE

5.1 Cleaning

WARNING: Electric Shock Hazard. Disconnect the equipment from the power supply before cleaning. Make sure that no liquid enters the interior of the instrument.

Attention: Do not use solvents, harsh chemicals, ammonia or abrasive cleaning agents.

The housing may be cleaned with a cloth dampened with a mild detergent if necessary.

5.2 Service Information

For troubleshooting common problems, download the user manual from www.Ohaus.com. If the issue persists, please visit our website for available support and service centers at <http://www.ohaus.com/en-us/support/service-centers>.

6. TECHNICAL DATA

Equipment Ratings:

- Indoor use only
- Altitude: 2000m
- Operating temperature: 0 to 40°C
- Humidity: Maximum relative humidity 80% for temperatures up to 31 °C decreasing linearly to 50% relative humidity at 40°C.
- Electrical supply: 12VDC, 0.5A. For use with certified or approved power supply, which must have a SELV and limited energy circuit output.
- Voltage fluctuations: Mains supply voltage fluctuations up to ±10% of the nominal voltage.
- Overvoltage category (Installation category): II
- Pollution degree: 2

Detailed product specifications can be found in the user manual and can be downloaded from www.Ohaus.com/products.

LIMITED WARRANTY

Ohaus products are warranted against defects in materials and workmanship from the date of delivery through the duration of the warranty period. During the warranty period Ohaus will repair, or replace any component(s) that proves to be defective at no charge, provided that the product is returned, freight prepaid, to Ohaus.

This warranty does not apply if the product has been damaged by accident or misuse, exposed to radioactive or corrosive materials, has foreign material penetrating to the inside of the product, or as a result of service or modification by other than Ohaus. In lieu of a properly returned warranty registration card, the warranty period shall begin on the date of shipment to the authorized dealer. No other express or implied warranty is given by Ohaus Corporation. Ohaus Corporation shall not be liable for any consequential damages.

As warranty legislation differs from state to state and country to country, please contact Ohaus or your local Ohaus dealer for further details.

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